

Human IL-1 β ELISA development kit

Product Code: 3415-1H-20

CONTENTS, development kit for 20 plates:

Vial 1 (green top)

Monoclonal antibody IL1 β -I (500 μ l)

Concentration: 1 mg/ml

Vial 2 (blue top)

Biotinylated monoclonal antibody IL1 β -II (250 μ l)

Concentration: 1 mg/ml

Vial 3 (white top)

Streptavidin-Horseradish Peroxidase (250 μ l)

Vial 4

Recombinant human IL-1 β standard (1 μ g)

To ensure total recovery of stated quantity, vials have been overfilled.

STORAGE:

Shipped at ambient temperature. On arrival box 1 should be stored refrigerated at 4-8°C and box 2 should be stored frozen at -20°C.

General

Intended use: For quantitative determination of native and recombinant human IL-1 β in solution, e.g. cell culture supernatant.

Serum/plasma samples: Please note that cytokine determinations in serum/plasma require the use of ELISA diluent (product code: 3652-D2) for dilution of samples, standard and detection antibody. The diluent prevents false positive read-outs which may be caused by interference of heterophilic antibodies found in plasma and serum. The ELISA diluent has been validated using serum/plasma from normal healthy human blood donors. Please note that heterophilic antibody interference in samples from human subjects with various diseases or other conditions has not been assessed. Please contact Mabtech for further information.

Reagents: Antibodies are supplied in sterile-filtered (0.2 μ m) PBS with sodium azide (0.02%). Streptavidin-HRP is supplied in PBS with 1% BSA and 0.002% Kathon CG.


Standard range: 3-300 pg/ml

Limit of detection: 1.5 pg/ml

Intra-assay variation: < 4%

Standard calibration: 1 ng of supplied standard equals 59 U of 86/680 NIBSC*-standard according to repeated calibrations. Calibration is batch-specific.

*National Institute of Biological Standards and Control, UK

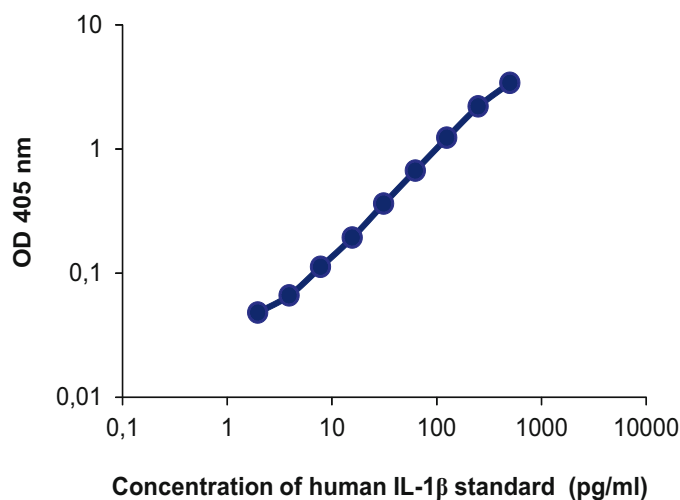

Guidelines for Human IL-1 β ELISA

- Day 1**
1. Coat a high protein binding ELISA plate with mAb IL1 β -I, diluted to 2 μ g/ml in PBS, pH 7.4, by adding 100 μ l/well. Incubate overnight at 4-8°C.
- Day 2**
2. Wash twice with PBS (200 μ l/well).
 3. Block plate by adding 200 μ l/well of PBS with 0.05% Tween 20 containing 0.1% BSA (incubation buffer). Incubate for 1 hour at room temperature.
 4. Wash five times with PBS containing 0.05% Tween20
 5. Prepare hIL-1 β standard by reconstituting contents of vial 4 in 1 ml PBS to a concentration of 1 μ g/ml. Leave at room temperature for 15 minutes and then vortex the tube. The stock solution should be used immediately or stored in aliquots at -20°C for future use. We recommend the aliquots not be refrozen after initial use. For the test, prepare dilutions of the stock using the standard range as a guideline.
 6. Add 100 μ l/well of samples or standards diluted in incubation buffer or ELISA diluent for serum/plasma samples and incubate for 2 hours at room temperature.
 7. Wash as in step 4.
 8. Add 100 μ l/well of mAb IL1 β -II-biotin at 1 μ g/ml in incubation buffer or ELISA diluent for serum/plasma samples. Incubate for 1 hour at room temperature.
 9. Wash as in step 4.
 10. Add 100 μ l/well of Streptavidin-HRP diluted 1:1000 in incubation buffer. Incubate for 1 hour at room temperature. **Please note that sodium azide used in buffers will inhibit HRP activity.**
 11. Wash as in step 4.
 12. Add 100 μ l/well of appropriate substrate solution
 13. Measure the optical density in an ELISA reader after suitable developing time. If required stop the reaction first.

MABTECH AB
Augustendalsvägen 19
Box 1233, SE-131 28 Nacka Strand
Sweden
Tel: +46 8 716 27 00
Fax: +46 8 716 27 01
E-mail: mabtech@mabtech.com

MABTECH Inc
M.E.B. 220, 3814 West Street
Cincinnati, OH 45227
USA
Toll free: +1 866 ELI-SPOT
Tel: +1 513 871 4500
Fax: +1 513 871 7353
E-mail: mabtech.usa@mabtech.com

MABTECH AUSTRALIA Pty Ltd
Australia
Tel: +61 3 9470 4704
Fax: +61 3 8678 3216
E-mail: mabtech.au@mabtech.com

MABTECH AB Büro Deutschland
Germany
Tel: +49 40 4135 7935
Fax: +49 40 4135 7945
E-mail: mabtech.de@mabtech.com

MABTECH AB Bureau de liaison
France
Tel: +33 (0)4 92 38 80 70
Fax: +33 (0)4 92 38 80 71
E-mail: mabtech.fr@mabtech.com

Updated on 2015-02-12

MABTECH

NOTE; for research use only.

MABTECH shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.



Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the above standards.

For more information about Mabtech services and our products, visit

www.mabtech.com