

Product Datasheet

PepPool: SARS-CoV-2 (SNMO), human

Product code:	3622-1
Contents:	The SARS-CoV-2 SNMO defined peptide pool contains 47 peptides from the human SARS-CoV-2 virus. The peptides are derived from the spike (S), nucleoprotein (N), membrane protein (M) and open reading frame (ORF)-3a, and ORF-7a proteins (O). The purity of the synthetic peptides range from 60-99%.
Applications:	The peptide pool is recommended for enumeration of cytokine secreting T cells specific for SARS-CoV-2 S, N, M, and O proteins with ELISpot/FluoroSpot. The peptide pool has been validated using human PBMC from COVID-19 convalescent individuals previously PCR-confirmed as SARS-CoV-2 positive.
Instructions:	<p>Sterile handling is recommended. Dissolve the lyophilized peptide pool by addition of 40 µl DMSO to the vial. Then add 85 µl PBS, mix and aliquote and store at -20°C or below. This stock solution will have a concentration of 200 µg/ml of each peptide.</p> <p>Dilute the stock solution 1:100 in cell culture medium to obtain 2 µg/ml of each peptide in the cell culture. Use the peptide pool in ELISpot and FluoroSpot assay for stimulation of 250,000-500,000 cells per well. Use the diluted peptide solution fresh.</p>
Storage:	Shipped at ambient temperature. Store frozen at -20°C or below upon receipt. After reconstitution, store aliquotes at -20°C or below. We recommend the aliquots not be refrozen after initial use.
Quantity:	One vial with 25 ug of each peptide.
	Reference: Peng Y, <i>et al.</i> Broad and strong memory CD4+ and CD8+ T cells induced by SARS-CoV-2 in UK convalescent COVID-19 patients. Nature Immunology, vol 21, Nov 2020.



Note; for research use only.

Mabtech shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.

Mabtech AB (Head Office)
Sweden
Tel: +46 8 716 27 00
mabtech@mabtech.com

Mabtech, Inc.
USA
Tel: +1 513 871-4500
mabtech.usa@mabtech.com

Updated on 2022-05-30



Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the standards ISO 9001:2015 & ISO 13485:2016.

www.mabtech.com

Peptides included in the SARS-CoV-2 SNMO defined peptide pool

Peptide	ID	Sequence	Epitope source
1	S-34	CTFEYVSQPFLMDLE	Spike
2	S-39	EFVFKNIDGYFKIYS	Spike
3	S-42	KHTPINLVRDLPOGF	Spike
4	S-43	NLVRDLPOGFSALEP	Spike
5	S-71	YAWNRRKRISNCVADY	Spike
6	S-77	GVSPTKLNLCFTNV	Spike
7	S-90	GGNYNYLYRFRKSN	Spike
8	S-91	YLYRFRKSNLKPFE	Spike
9	S-103	VVLSFELLHAPATVC	Spike
10	S-106	GPKKSTNLVKNKCVN	Spike
11	S-145	SVTTEILPVSMTKTS	Spike
12	S-150	STECNLLLQYGSFC	Spike
13	S-151	NLLLQYGSFCTQLNR	Spike
14	S-161	NFSQILPDPSPSKR	Spike
15	S-174	TDEMIAQYTSALLAG	Spike
16	S-235	GINASVVNIQKEIDR	Spike
17	S-240	LIDLQELGKYEQYI	Spike
18	S-242	YEQYIKWPWYIWLGF	Spike
19	NP-1	MSDNGPQNQRNAPRITF	Nucleoprotein
20	NP-2	NQRNAPRITFGGPSDSTG	Nucleoprotein
21	NP-12	DQIGYYRRATRRIR	Nucleoprotein
22	NP-15	MKDLSRWYFYFL	Nucleoprotein
23	NP-16	LSPRWYFYFLGTGPEAGL	Nucleoprotein
24	NP-46	AFFGMSRIGMEVTPSGTW	Nucleoprotein
25	NP-47	GMEVTPSGTWLTYTGAIK	Nucleoprotein
26	NP-48	TWLTYTGAIKLDDKPNF	Nucleoprotein
27	NP-50	PNFKDQVILLNKHIDAYK	Nucleoprotein
28	NP-51	LLNKHIDAYKTFPPEPK	Nucleoprotein
29	M-19	LLESELVIGAVILRGHLR	Membrane protein
30	M-20	GAVILRGHLRIAGHHLGR	Membrane protein
31	M-21	LRIAGHHLGRCDIKDLPK	Membrane protein
32	M-23	PKEITVATSRTLSYYKL	Membrane protein
33	M-24	TSRTLSYYKLGASQRVA	Membrane protein
34	M-28	IGNYKLNTDHSSSDNIA	Membrane protein
35	ORF-3a-20	YFLCWHTNCYDYCIPY	ORF 3a
36	ORF-3a-27	KDCVVLSYFTSDYYQLY	ORF 3a
37	ORF-3a-28	YFTSDYYQLYSTQLSTDTGV	ORF 3a
38	ORF-3a-30	GVEHVTFYFNKIVDEPEEH	ORF 3a
39	ORF-7a-2	LITLATCELYHYQECVR	ORF 7a
40	ORF-7a-7	FHPLADNKFALTCFSTQF	ORF 7a
41	ORF-7a-10	DGVKHVYQLRARSVSPKL	ORF 7a
42	N-E-01	ILLNKHID	Nucleoprotein
43	N-E-03	MEVTPSGTWL	Nucleoprotein
44	S-E-19	QLIRAAEIRASANLAATK	Spike
45	N	SPRWYFYFL	Nucleoprotein
46	N	YLGTGPEAGL	Nucleoprotein
47	N	YYLGTGPEA	Nucleoprotein