

Product Datasheet

FluoroSpot enhancer

Product code:	3641-F10
Contents:	Ready-to-use solution with 0.002% Kathon CG as preservative.
Applications:	The Fluorescence enhancer will increase spot quality in FluoroSpot assays. Please note that the Fluorescence enhancer is optimized for use with FluoroSpot assays from Mabtech.
Instructions:	The Fluorescence enhancer should be added to the plate after removal of the fluorophore conjugate and a 5 x PBS wash. Add 50 µl per well of Fluorescence enhancer and leave for 15 minutes at room temperature. Empty and remove residual enhancer by firmly tapping the plate against clean paper towels.
Storage:	Shipped at ambient temperature. Store at 4-8°C upon receipt. 120 ml



Note; for research use only.

Mabtech shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.

Mabtech AB (Head Office)
Sweden
Tel: +46 8 716 27 00
mabtech@mabtech.com

Mabtech, Inc.
USA
Tel: +1 513 871-4500
mabtech.usa@mabtech.com

Updated on 2022-05-31



Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the standards ISO 9001:2015 & ISO 13485:2016.

www.mabtech.com