

Product Datasheet

Apo ELISA buffer

Product code:	3652-M2
Contents:	Protein-containing buffer, 5x concentrated solution, with 0.002% Kathon CG as preservative.
Applications:	To prevent interference by heterophilic antibodies in serum and plasma samples in ELISA. Recommended for use with apolipoprotein antibodies and ELISA kits from Mabtech. Please note that the Apo ELISA buffer is optimized for use with antibodies from Mabtech.
Instructions:	<p>The Apo ELISA buffer is supplied as a 5x concentrated solution. Before use, dilute five times with distilled or deionized water. The Apo ELISA buffer should be used for dilution of samples, ELISA standard and detection antibody. All samples should be diluted at least 2x. Do not use undiluted serum/plasma samples in the assay. Sample dilution can be made either in separate tubes or directly in the plate. If sample dilution is made directly in the plate, Apo ELISA buffer should be added prior to addition of the sample.</p> <p>Important! Do not use the Apo ELISA buffer for dilution of Streptavidin conjugates. We recommend the use of an incubation buffer e.g. PBS with 0.05% Tween 20 containing 0.1% BSA for dilution of conjugates.</p>
Storage:	Shipped at ambient temperature. Store at 4-8°C upon receipt. Adjust to room temperature before use.



Note; for research use only.

Mabtech shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.

Mabtech AB (Head Office)
Sweden
Tel: +46 8 716 27 00
mabtech@mabtech.com

Mabtech, Inc.
USA
Tel: +1 513 871-4500
mabtech.usa@mabtech.com

Updated on 2022-05-31



Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the standards ISO 9001:2015 & ISO 13485:2016.

www.mabtech.com