

Human IgM ELISA development kit

Product Code: 3840-1AD-6

CONTENTS:

Vial 1 (red top)

Monoclonal anti-IgM antibody (300 µl)

Concentration: 0.5 mg/ml

Vial 2 (yellow top)

ALP-conjugated anti-IgM antibody (80 µl)

Vial 3

Lyophilised human IgM standard

To ensure total recovery of stated quantity, vials have been overfilled.

STORAGE:

Shipped at ambient temperature. On arrival box 1 should be stored refrigerated at 4-8°C and box 2 should be stored frozen at -20°C.

General


Intended use: For quantitative determination of native human IgM in serum and plasma.

Reagents: Anti-IgM monoclonal antibody is supplied in sterile-filtered (0.2 μm) PBS with sodium azide (0.02%). ALP-conjugated anti-IgM antibody is supplied in 0.1 M Tris-buffer with 0.002% Kathon CG.

Recommended standard dilution: 0.1-500 ng/ml

Standard calibration: 1 μg of supplied standard equals 233 mU NIBSC* standard. Calibration is batch-specific.

* National Institute of Biological Standards and Control, UK.

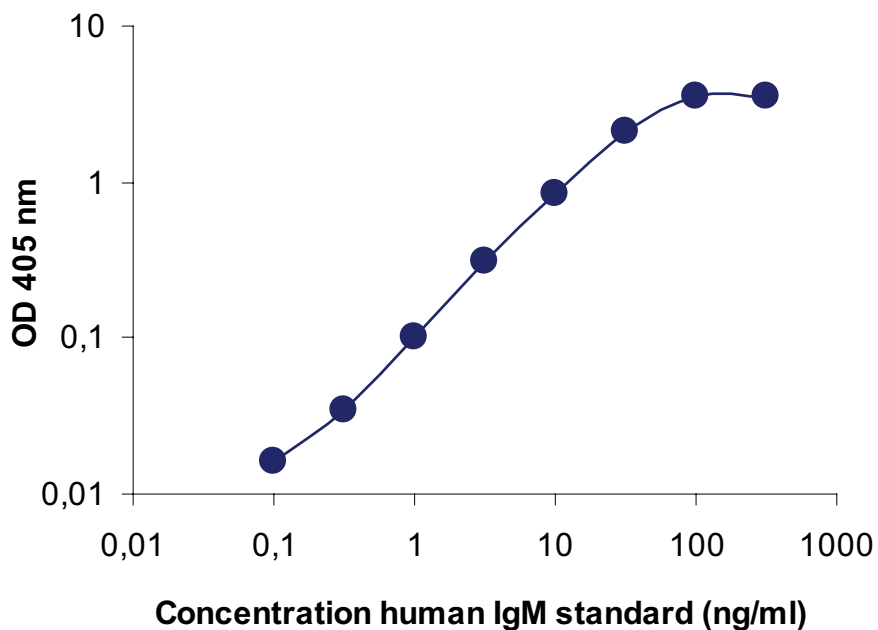

Guidelines for Human IgM ELISA

- Day 1**
1. Coat a high protein binding ELISA plate with anti-IgM monoclonal antibody, diluted to 2 µg/ml in PBS, pH 7.4, by adding 100 µl/well. Incubate overnight at 4-8°C.
- Day 2**
2. Wash twice with PBS (200 µl/well).
 3. Block plate by adding 200 µl/well of PBS with 0.05% Tween 20 (PBS-Tween) containing 0.1% BSA (incubation buffer*). Incubate for 1 hour at room temperature.
 4. Wash five times with PBS-Tween.
 5. Prepare human IgM standard by reconstituting contents of vial 3 in 500 µl PBS to make up a stock solution of 50 µg/ml. The stock solution should be used immediately or stored in aliquots at -20°C for future use. We recommend the aliquots not be refrozen after initial use. For the test, prepare dilutions of the stock using the standard range as a guideline.
 6. Add 100 µl/well of samples or standards diluted in incubation buffer. It is recommended to make serial dilutions of samples to obtain results within the standard range. Incubate for 2 hours at room temperature.
 7. Wash as in step 4.
 8. Add 100 µl/well of anti-IgM-ALP diluted 1:1000 in incubation buffer. Incubate for 1 hour at room temperature.
 9. Wash as in step 4.
 10. Add 100 µl/well of appropriate substrate solution e.g. p-nitrophenyl-phosphate (pNPP).
 11. Measure the optical density (405 nm for pNPP) in an ELISA reader after suitable developing time.

* The same buffer is used for blocking and for dilution.

MABTECH AB
Augustendalsvägen 19
Box 1233, SE-131 28 Nacka Strand
Sweden
Tel: +46 8 716 27 00
Fax: +46 8 716 27 01
E-mail: mabtech@mabtech.com

MABTECH Inc
M.E.B. 220, 3814 West Street
Cincinnati, OH 45227
USA
Toll free: +1 866 ELI-SPOT
Tel: +1 513 871 4500
Fax: +1 513 871 7353
E-mail: mabtech.usa@mabtech.com

MABTECH AUSTRALIA Pty Ltd
Australia
Tel: +61 3 9470 4704
Fax: +61 3 8678 3216
E-mail: mabtech.au@mabtech.com

MABTECH AB Büro Deutschland
Germany
Tel: +49 40 4135 7935
Fax: +49 40 4135 7945
E-mail: mabtech.de@mabtech.com

MABTECH AB Bureau de liaison
France
Tel: +33 (0)4 92 38 80 70
Fax: +33 (0)4 92 38 80 71
E-mail: mabtech.fr@mabtech.com

MABTECH

NOTE; for research use only.

MABTECH shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.



Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the above standards.

For more information about Mabtech services and our products, visit

www.mabtech.com