Human IgM ELISA development kit

Product Code: 3880-1AD-6

CONTENTS: development kit for 6 plates

Vial 1 (red top)

Monoclonal antibodies MT11/12 (300 μl)

Concentration: 0.5 mg/ml

Vial 2 (yellow top)

Monoclonal antibody MT22-ALP (80 µl)

Vial 3

Lyophilised human IgM standard

To ensure total recovery of stated quantity, vials have been overfilled.

STORAGE:

Shipped at ambient temperature. On arrival box 1 should be stored refrigerated at 4-8°C and box 2 should be stored frozen at -20°C.

General


Intended use: For quantitative determination of native human IgM in serum, plasma or cell supernatant.

Reagents: MAbs MT11/12 are supplied in sterile-filtered (0.2 μ m) PBS with sodium azide (0.02%). MT22-ALP is supplied in 0.1 M Tris-buffer with 1% BSA and 0.002% Kathon.

Standard range: 0.2-200 ng/ml

Standard calibration: 1 µg of supplied standard equals 233 mU NIBSC* standard. Calibration is batch-specific.

* National Institute of Biological Standards and Control, UK.

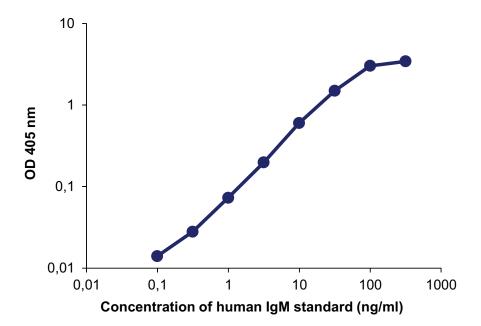

Guidelines for Human IgM ELISA

- Day 1 1. Coat a high protein binding ELISA plate with mAbs MT11/12, diluted to 2 μg/ml in PBS, pH 7.4, by adding 100 μl/well. Incubate overnight at 4-8°C.
- **Day 2** 2. Wash twice with PBS (200 μl/well).
 - 3. Block plate by adding 200 µl/well of PBS with 0.05% Tween 20 (PBS-Tween) containing 0.1% BSA (incubation buffer*). Incubate for 1 hour at room temperature.
 - 4. Wash five times with PBS-Tween.
 - 5. Prepare human IgM standard by reconstituting contents of vial 3 in 500 µl PBS to make up a stock solution of 50 µg/ml. The stock solution should be used immediately or stored in aliquots at -20°C for future use. We recommend the aliquots not be refrozen after initial use. For the test, prepare dilutions of the stock using the standard range as a guideline.
 - 6. Add 100 µl/well of samples or standards diluted in incubation buffer. It is recommended to make serial dilutions of samples to obtain results within the standard range. Incubate for 2 hours at room temperature.
 - 7. Wash as in step 4.
 - 8. Add 100 µl/well of MT22-ALP diluted 1:1000 in incubation buffer. Incubate for 1 hour at room temperature.
 - 9. Wash as in step 4.
 - 10. Add 100 μl/well of appropriate substrate solution e.g. p-nitrophenyl-phosphate (pNPP).
 - 11. Measure the optical density (405 nm for pNPP) in an ELISA reader after suitable developing time.
 - * The same buffer is used for blocking and for dilution.



Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the standards ISO 9001:2015 & ISO 13485:2016.





The products are for research use only.

MABTECH shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages there from.

Mabtech AB (Head Office)

Sweden Tel: +46 8 716 27 00 mabtech@mabtech.com Mabtech, Inc.

USA Tel: +1 513 871-4500 mabtech.usa@mabtech.com